Metallpyknometer

Metal pyknometer

ZPM 3030

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The density is an essential quality characteristic, for example it allows recognising errors regarding the composition and gives indications of the purity of a product.

The test principle is based on the weighing of a given volume of the liquid under test, as it is defined by the pyknometer. From the difference of the masses of the filled and the empty pyknometer as well as the known filling volume the density is calculated.

Application/application areas

- laboratory apparatus for the quality control and for experimental purposes
- for very different branches of industry, e.g. for the paint, cosmetics, food, cleaning agents industries and the chemical industry in general, as well as for technical high schools, universities and testing institutes

Standard extent of delivery

- 1 Pyknometer, consisting of a cylindrical beaker with a lid having a hole
- 1 certificate of manufacturer

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Handling

- weigh the carefully cleaned empty pyknometer plus lid and record the mass
- bring the pyknometer and the substance under test to 23 °C ± 0,5 °C, unless no other test temperature is specified or agreed
- fill the pyknometer with the substance under test and place the lid firmly in position without tilting, whilst avoiding inclusion of air bubbles
- remove any excess liquid from the lid using a clean and absorbent cloth
- weigh the filled pyknometer
- calculate the density, using the following formula:
density = mass (g) of filled pyknometer minus mass (g) of empty pyknometer, divided by the volume (ml) of the pyknometer
- do not use sharp-edged objects when cleaning the pyknometer in order to avoid damaging, but a suitable solvent

Technical specification

	filling volume	height/diameter	weight
ZPM 3030.50	50 ml	35 mm x Ø 54 mm	144.1 g
ZPM 3030.100	100 ml	63 mm x Ø 54 mm	174.4 g

material: stainless steel
 manufacturing: ± 0,2 %, related to the volume;
 tolerance: as option with certificate of calibration ± 0,1 %
 standards: ASTM D 1475, EN ISO 2811-1,
 SNV 37100, VDA 621-103
 warranty: 2 years

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ZPM 3030 Pyknometer distributed by
 Zehntner GmbH Testing Instruments.
 Subject to technical changes.

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